

## Your distributor

### Europe

#### Austria

VWR International GmbH  
Graumanngasse 7  
1150 Wien  
Tel.: 01 97 002 0  
Fax: 01 97 002 600  
E-mail: info@at.vwr.com

#### Belgium

VWR International bvba  
Researchpark Haasrode  
2020  
Geldenaaksebaan 464  
3001 Leuven  
Tel.: 016 385 1011  
Fax: 016 385 385  
E-mail:  
customerservice@be.vwr.com

#### Denmark

VWR - Bie & Berntsen  
Transformervej 8  
2730 Herlev  
Tel.: 43 86 87 88  
Fax: 43 86 87 90  
E-mail: info@dk.vwr.com

#### Finland

VWR International Oy  
Valimotie 9  
00380 Helsinki  
Tel.: 09 80 45 51  
Fax: 09 80 45 52 00  
E-mail: info@fi.vwr.com

#### France

VWR International S.A.S.  
Le Périgares – Bâtiment B  
201, rue Carnot  
94126 Fontenay-sous-Bois  
cedex  
Tel.: 0 825 02 30 30 (0.15  
EUR TTC/mn)  
Fax: 0 825 02 30 35 (0.15  
EUR TTC/mn)  
E-mail: info@fr.vwr.com

#### Germany

VWR International GmbH  
Hilpertstrasse 20a  
D - 64295 Darmstadt  
Tel.: 0180 570 20 00\*  
Fax: 0180 570 22 22\*  
E-mail: info@de.vwr.com  
\*0.14 €/Min. aus d. dt. Festnetz,  
Mobilfunk max. 0.42 €/Min.

#### Hungary

VWR International Kft.  
Simon László u. 4.  
4034 Debrecen  
Tel.: (52) 521-130  
Fax: (52) 470-069  
E-mail: info@hu.vwr.com

#### Ireland / Northern Ireland

VWR International Ltd /  
VWR International (North-  
ern Ireland) Ltd  
Orion Business Campus  
Northwest Business Park  
Ballycoolin  
Dublin 15  
Tel.: 01 88 22 222  
Fax: 01 88 22 333  
E-mail: sales@ie.vwr.com

#### Italy

VWR International PBI  
S.r.l.  
Via San Giusto 85  
20163 Milano (MI)  
Tel.: 02-3320311/02-  
487791  
Fax: 800 152999/02-  
40090010  
E-mail: info@it.vwr.com  
info@internationalpbi.it

#### The Netherlands

VWR International B.V.  
Postbus 8198  
1005 AD Amsterdam  
Tel.: 020 4808 400  
Fax: 020 4808 480  
E-mail: info@nl.vwr.com

#### Norway

VWR International AS  
Haavard Martinsens vei 30  
0978 Oslo  
Tel.: 02290  
Fax: 815 00 940  
E-mail: info@no.vwr.com

#### Poland

Labart Sp. z o.o.  
A VWR International Com-  
pany  
Limbowa 5  
80-175 Gdansk  
Tel.: 058 32 38 200 do 204  
Fax: 058 32 38 205  
E-mail: labart@pl.vwr.com

#### Portugal

VWR International - Mate-  
rial de Laboratório, Lda  
Edifício Neopark  
Av. Tomás Ribeiro, 43- 3  
□  
2790-221 Carnaxide  
Tel.: 21 3600 770  
Fax: 21 3600 798/9  
E-mail: info@pt.vwr.com

#### Spain

VWR International Eurolab  
S.L.  
C/ Tecnología 5-17  
A-7 Llinars Park  
08450 - Llinars del Vallès  
Barcelona  
Tel.: 902 222 897  
Fax: 902 430 657  
E-mail: info@es.vwr.com

#### Sweden

VWR International AB  
Fagerstagatan 18a  
163 94 Stockholm  
Tel.: 08 621 34 00  
Fax: 08 621 34 66  
E-mail: info@se.vwr.com

#### Switzerland

VWR International AG  
Lerzenstrasse 16/18  
8953 Dietikon  
Tel.: 044 745 13 13  
Fax: 044 745 13 10  
E-mail: info@ch.vwr.com

#### UK

VWR International Ltd  
Customer Service Centre  
Hunter Boulevard  
Magna Park  
Lutterworth  
Leicestershire  
LE17 4XN  
Tel.: 0800 22 33 44  
Fax: 01455 55 85 86  
E-mail: uk-  
sales@uk.vwr.com

### Asia Pacific

#### China

VWR International China  
Co., Ltd  
Suite 3B02, Qilai Building,  
No. 889  
Yishan Road  
Shanghai 200233, China  
Tel.: +86-21 521 388 22  
Fax: +86-21 521 33 933  
E-mail: sa-  
les\_china@vwr.com

#### India

VWR Lab Products Pvt.  
Ltd  
2nd Floor, Front Wing,  
135/12, Brigade Towers  
Brigade Road  
Bangalore 560025 India  
Tel.: +91-2522-  
647911/922  
(Mumbai)  
Tel.: +91-80-41117125/26  
(Bangalore)  
Fax: +91-80-41117120  
E-mail:  
vwr\_india@vwr.com

#### Singapore

VWR Singapore Pte Ltd  
18 Gul Drive  
Singapore 629468  
Tel.: +65 6505 0760  
Fax: +65 6264 3780  
E-mail: sales@sg.vwr.com

H35 0.01 SC-07 VWR

as of version V2.9

VWR

EN

Quick-Response Double / Difference Thermometer  
For Thermocouple Sensor Type J, K, N, S, T

Operating Manual

TD 241

European Catalogue Number(s):  
Euro Plug: 620-1935

Issued: 01.04.2013

Go to [vwr.com](http://vwr.com) for the latest news, special offers and details of  
your local VWR distributor

CE

Made in  
Germany

WEEE-Reg.-Nr. DE93889386

## Index

|      |  |    |
|------|--|----|
| 1    | DESIGNATED USE                                       | 2  |
| 2    | GENERAL NOTE   | 2  |
| 3    | OPERATING AND MAINTENANCE ADVICE                     | 3  |
| 4    | SAFETY REQUIREMENTS                                  | 3  |
| 5    | DISPLAY AND CONTROL ELEMENTS                         | 4  |
| 5.1  | DISPLAY ELEMENTS                                     | 4  |
| 5.2  | PUSHBUTTONS  | 5  |
| 5.3  | CONNECTIONS  | 5  |
| 6    | CONFIGURATION  | 5  |
| 7    | MEASURING TEMPERATURE WITH THERMOELEMENTS            | 7  |
| 8    | NOTES TO SPECIAL FUNCTIONS                           | 7  |
| 8.1  | DISPLAY RESOLUTION ('RESOLUTION')                    | 7  |
| 8.2  | ZERO DISPLACEMENT ('OFFSET')                         | 7  |
| 8.3  | DISPLAY CORRECTION FACTOR ('CORR')                   | 7  |
| 8.4  | TARE FUNCTION  | 7  |
| 8.5  | BASE ADDRESS ('ADR.')                                | 7  |
| 9    | ERROR AND SYSTEM MESSAGES                            | 8  |
| 10   | THE SERIAL INTERFACE                                 | 9  |
| 10.1 | THE FOLLOWING INTERFACE FUNCTIONS WILL BE SUPPORTED: | 9  |
| 11   | SPECIFICATION  | 10 |
| 12   | TECHNICAL SERVICE                                    | 11 |
| 13   | WARRANTY   | 11 |
| 14   | EQUIPMENT DISPOSAL                                   | 11 |

## 1 Designated Use

This device is designed for temperature measurement.

You can plug in up to 2 thermocouple sensors with Minitur-DIN-plug.

If you use two sensors, you can measure two different temperatures as well as the difference between them.

The device features a lot of useful features as for example min-/max-alarm, hold, correction factor for surface measurement, real-time clock, interface etc.

## 2 General Note

Read this document carefully and get used to the operation of the device before you use it. Keep this document within reach for consulting in case of doubt.

## 3 Operating and Maintenance Advice

- When to replace battery:  
If  $\Delta$  and 'bAt' are shown in the lower display the battery has been used up and needs to be replaced. The device will, however, operate correctly for a certain time.  
If 'bAt' is shown in the upper display the voltage is too low to operate the device; the battery has been completely used up.  
*Please note: The battery has to be taken out, when storing device above 50°C.  
We recommend to take out battery if device is not used for a longer period of time.*
- Treat device and sensor carefully. Use only in accordance with above specification. (do not throw, hit against etc.). Protect plug and socket from soiling.
- To disconnect sensor thermoelement plug do not pull at the cable but at the plug.  
If plug is entered incorrectly the connecting pins of the plug can be damaged. => Plug can no longer be used and connecting cable needs to be replaced.
- Selection of types of thermoelements:  
Prior to carrying out a measurement make sure to check if device is set to the thermoelement type used (type is shown on the display shortly after device has been switched on). Unless the correct thermoelement is set, temperature measurements will be incorrect.
- Mains operation:  
When using a power supply device please note that operating voltage has to be 10.5 to 12 V DC.  
Do not apply overvoltage!! Cheap 12V-power supply devices often have excessive no-load voltage. We, therefore, recommend using regulated voltage power supply devices. Trouble-free operation is guaranteed by our power supply devices. Trouble-free operation is guaranteed by our power supply, PS 100.  
Prior to connecting the plug power supply device with the mains supply make sure that the operating voltage stated at the power supply device is identical to the mains voltage.

## 4 Safety Requirements

This device has been designed and tested in accordance with the safety regulations for electronic devices. However, its trouble-free operation and reliability cannot be guaranteed unless the standard safety measures and special safety advises given in this manual will be adhered to when using the device.

- Trouble-free operation and reliability of the device can only be guaranteed if the device is not subjected to any other climatic conditions than those stated under "Specification".
- If the device is transported from a cold to a warm environment condensation may cause in a failure of the function. In such a case make sure the device temperature has adjusted to the ambient temperature before trying a new start-up.
- If device is to be connected to other devices (e.g. via serial interface) the circuitry has to be designed most carefully. Internal connection in third party devices (e.g. connection GND and earth) may result in not-permissible voltages impairing or destroying the device or another device connected.  
**Warning:** If device is operated with a defective mains power supply (short circuit from mains voltage to output voltage) this may result in hazardous voltages at the device (e.g. sensor socket, serial interface).
- If there is a risk whatsoever involved in running it, the device has to be switched off immediately and to be marked accordingly to avoid re-starting.  
Operator safety may be a risk if:
  - there is visible damage to the device
  - the device is not working as specified
  - the device has been stored under unsuitable conditions for a longer time.
 In case of doubt, please return device to manufacturer for repair or maintenance.
- Warning:** Do not use these products as safety or emergency stop devices or in any other application where failure of the product could result in personal injury or material damage.  
Failure to comply with these instructions could result in death or serious injury and material damage.

## 5 Display and control elements

### 5.1 Display elements

#### Display with two sensors connected:



① **Main display:** possible views:

- CH1: sensor 1
- CH2: sensor 2
- DIF: sensor 1 - sensor 2

② **Secondary display:** possible views:

- CH1: sensor 1
- CH2: sensor 2
- DIF: sensor 1 - sensor 2

The desired view can be selected by pressing the -key.

#### Display with one sensor connected:



① **Main display:** possible views:

- CH1: sensor 1
- CH2: sensor 2

Device automatically detects to which socket sensor is connected.

#### Special display elements:



① **Min/Max/Hold:** shows if a min., max. or hold value is displayed in either the main or the secondary display.

② **Offset arrow:** indicates that zero point offset (offset) is activated

③ **Corr arrow:** indicates that correction factor is activated

④ **Warning triangle:** indicates a low battery, full logger storage, etc.

⑤ **Tare arrow:** indicates that tare function is activated (for 'DIF'-display when working with two sensors only)

#### Display at reboot:

After switching on the instrument performs a segment test of 2 seconds. Afterwards some configurations are displayed sequentially: thermocouple type and if activated: offset and scale settings. (p.r.t Chapter 6)

*Tip:* The display of these settings can be aborted by pressing a key after the segment test (keys 2 - 6).

### 5.2 Pushbuttons



**On/off key**

**min/max when taking measurements:**

press shortly: min. or max. measuring value will be displayed

press for 1 sec.: the min. or max. value will be deleted up/down for config. or selection for logger operation: to enter values, or change settings

**Tare:** (for 'DIF'-display when working with two sensors only)

press shortly: the difference between CH1-CH2 will be set to zero.

press for 1 sec.: tare function will be deactivated

**Set/Menu:**

press (Set) shortly in case of 2 sensors: the values displayed (CH1, CH2, DIF) can be selected  
press (Menu) for 2 sec.: configuration will be activated

**Store/Quit:**

Measurement: Hold current measuring value ('HLD' in display)

Set/Menu: Acknowledge setting, return to measuring

### 5.3 Connections



The mains socket is located at the left side of the instrument.

## 6 Configuration

For configuration of the device press -key for 2 seconds: the main menu of the configuration will be called up. Use key to select a sub-menu, use the key to actually go into the selected sub-menu and to change parameters.

Use the keys and to set the individual value for the parameter. Press the key again to memorize the changes and to change to the main menu. Use key to leave the configuration.



#### 'Typ': Selection of Thermoelement Type\*1

- |  |        |                     |
|--|--------|---------------------|
|  | ni.cr: | Type K: NiCr-Ni     |
|  | n:     | Type N: NiCrSi-NiSi |
|  | S:     | Type S: Pt10Rh-Pt   |
|  | t:     | Type T: Cu-CuNi     |
|  | J:     | Type J: Fe-CuNi     |



#### 'Resolution': Selection of Display Resolution

- |  |       |                                      |
|--|-------|--------------------------------------|
|  | 1°:   | Resolution 1°C                       |
|  | 0.1°: | Resolution 0.1°C                     |
|  | Auto: | Resolution is selected automatically |



### 'Unit': Selection of Temperature Unit °C /°F\*1



°C: All temperature values in degrees Celsius



°F: All temperature values in degrees Fahrenheit



### 'Corr': Selection of Display Correction Factor \*1



0.950 ... 1.200:

The temperature value (referring to 0°C or 32°F) will be multiplied by this factor. This factor will be applied to both sensor values.



off:

Factor is deactivated (=1.000)



### 'Offset': Zero Displacement Channel 1 \*1



-10.0°C ... 10.0°C

The zero point of the measurement of channel 1 will be displaced by this value.



or  
-18.0°F ... 18.0°F

off:

Zero point displacement is deactivated (=0.0°)



### 'Offset': Zero Displacement Channel 2 \*1



-10.0°C ... 10.0°C

The zero point of the measurement of channel 2 will be displaced by this value.



or  
-18.0°F ... 18.0°F

off:

Zero point displacement is deactivated (=0.0°)



### 'Power.off': Selection of Power-Off Delay



1...120:

Power-off delay in minutes. Device will be automatically switched off as soon as this time has elapsed if no key is pressed/no interface communication takes place. (automatically deactivated for cyclic loggers)



off:

automatic power-off function deactivated (continuous operation, e.g. in case of mains operation)



### 'Adresse': Selection of Base Address'



01, 11, 21, ..., 91:

Base address for interface communication. Channel 1 will be addressed by the set base address, channel 2 and 3 will have the following addresses. (Example: base address 21 - channel 1 = 21, channel 2 = 22, channel 3 = 23)



Using the interface converter it is possible to connect several devices to a single interface. As a precondition the base addresses of all devices must not be identical. In case several devices are connected via one interface make sure to configure the base addresses accordingly.

## 7 Measuring Temperature With Thermoelements

- Make sure that the correct type of thermoelement is set (p.r.t. "Configuration of The Instrument")! Otherwise a wrong temperature will be displayed. The measuring device has been optimised for measurements with a type K probe.
- If other thermoelements but type K (NiCr-Ni) are used already the smallest temperature difference between plug and instrument will result in measuring errors. Therefore wait for the temperatures to adjust after plug-in or touching a sensor plug (depending on temperature ~15min.)
- Thermoelements are suitable to measure within a large temperature range. But keep in mind the allowed temperature range of your specific temperature probe!

## 8 Notes to Special Functions

### 8.1 Display Resolution ('Resolution')

Standard setting: 'Auto', i.e. the device automatically switches over to the optimum resolution between 1° and 0.1°. If temperatures to be measured are near the switching threshold, a fixed resolution may be better, e.g. for easy recording.

In such a case please select the optimum resolution manually.

### 8.2 Zero Displacement ('Offset')

A zero displacement can be carried out for each of the two channels CH1 or CH2.

**temperature displayed = temperature measured - offset**

Standard setting: 'off' = 0.0°, i.e. no zero displacement will be carried out. The zero displacement is mainly used to compensate for sensor deviations. Unless 'off' is set, this value will be displayed shortly after the device is switched on; during operation it will be identified by means of the offset arrow in the display.

### 8.3 Display Correction Factor ('Corr')

This factor is applied to both sensor channels.

**temperature displayed [°C] = temperature measured [°C] \* Corr**

**or temperature displayed [°F] = (temperature measured [°F]-32°F) \* Corr + 32°F**

Standard setting: 'off' = 1.000

This factor is used to compensate for losses of transfer in case of surface measurements, occurring if the object to be measured is extremely hot but will be cooled by lower ambient temperatures. The same can be true for sensors with a large mass. Unless 'off' is set, this value will be displayed shortly after the device is switched on; during operation it will be identified by means of the Corr-arrow in the display.

### 8.4 Tare Function




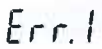

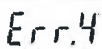

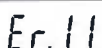
The tare function is used to set the DIF-display to zero, which is especially important when monitoring temperature differences. If the tare key is pressed for more than 2 seconds DIF = CH1-CH2 will be restored. This function as well as the DIF-display can only be used if two sensors are connected. If the tare function is activated, this will be shown by the tare arrow in the display.

### 8.5 Base Address ('Adr.')

Using the interface converter it is possible to connect several instruments to a single interface. As a precondition the base addresses of all devices must not be identical. In case several devices will be connected via one interface make sure to configure the base addresses accordingly. Channel 1 will be addressed by the base address set, channels 2 and 3 will have the following addresses.

(Example: base address 21 - channel 1 = 21, channel 2 = 22, channel 3 = 23)

## 9 Error And System Messages

| Error or system messages  | Description / Reason  | Remedy   |
|---|---|--|
|    | No sensor connected<br>sensor/cable damaged                                 | Connect one sensor at least<br>Replace sensor/cable  |
|    | Low battery voltage<br>device will only continue operation for a short time | replace battery  |
|    | Low battery voltage<br>- If mains operation: wrong voltage                  | replace battery<br>replace power supply, if fault continues to exist:<br>device damaged  |
| No display<br>or<br>mazy characters   | - Battery voltage too low   | replace battery  |
|   | - If mains op.: power supply defective or wrong voltage/polarity            | check/replace mains supply   |
|   | - System error  | disconnect battery or power supply, wait for a short time, re-connect  |
|    | - Device defective  | return to manufacturer for repair  |
|   | Values exceeding measuring range  | Are there any values exceeding the measuring range specified?<br>->temperature too high<br>Display resolution set to 0.1°?<br>-> set to 'Auto'   |
|    | Sensor/cable defective  | -> replace probe   |
|   | Values below measuring range  | Are there any values below the measuring range specified?<br>-> temperature too low<br>Display resolution set to 0,1°? -> set to 'Auto' for DIF-channel and difference<-2040°<br>-> temp. difference too low |
|   | Sensor/cable defective  | -> replace probe   |
|   | Values below display range  | for DIF-channel and difference <-1999°:<br>exchange sensors<br>CH1<->CH2 (watch out for offset setting!)   |
|  | System fault  | switch on again: if fault continues to exist, device is damaged<br>-> return to manufacturer for repair  |
|   | Instrument not within working temperature                                   | keep working temperature in between -25...50°C   |
|  | Value could not be calculated   | One measuring variable required for calculation is missing (no sensor) or incorrect (overflow/underflow)   |

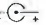
## 10 The serial interface

All measuring data and settings of the device can be read and changed by means of the serial interface and a suitable electrically isolated interface adapter.  
In order to avoid faulty transmission, we have designed elaborate security measures for interface communication.

### 10.1 The following interface functions will be supported:

| Channel |   |   | DII-Code | Name / function                |
|---------|---|---|----------|--------------------------------|
| 1       | 2 | 3 |          |                                |
| x       | x | x | 0        | Read nominal value             |
| x       | x | x | 3        | Read system status             |
| x       | x | x | 6        | Read min. value                |
| x       | x | x | 7        | Read max. value                |
| x       |   |   | 12       | Read ID-no.                    |
| x       |   |   | 174      | Delete min. value              |
| x       |   |   | 175      | Delete max. value              |
| x       |   |   | 194      | Set display unit               |
| x       |   |   | 195      | Set decimal point in display   |
| x       | x | x | 199      | Read meas. type in display     |
| x       | x | x | 200      | Read min. display range        |
| x       | x | x | 201      | Read max. display range        |
| x       | x | x | 202      | Read unit of display           |
| x       | x | x | 204      | Read decimal point of display  |
| x       |   |   | 208      | Read channel count             |
| x       | x |   | 216      | Read offset correction         |
| x       | x |   | 217      | Set offset correction          |
| x       |   |   | 218      | Read corr. factor (1000..1200) |
| x       |   |   | 219      | Set corr. factor (1000..1200)  |
| x       |   |   | 240      | Reset unit                     |
| x       |   |   | 254      | Read program identification    |

## 11 Specification

|                                     |   |   |
|-------------------------------------|---|---|
| <b>Thermoelements:</b>              | J, K, N, S, T   |   |
| <b>Resolution:</b>                  | 0.1°C or 1°C  | 0.1°F or 1°F                            |
| <b>Measuring range:</b>             |   |   |
| <b>Type K (NiCr-Ni):</b>            | -199.9 ... +999.9°C or -220 ... +1372°C   | -199.9 ... +999.9°F or -364 ... +2500°F |
| <b>Type J (Fe-CuNi):</b>            | -120.0 ... +700.0°C or -200 ... +1100°C   | -184.0 ... +999.9°F or -328 ... +2012°F |
| <b>Type N (NiCrSi-NiSi):</b>        | -199.9 ... +999.9°C or -200 ... +1300°C   | -199.9 ... +999.9°F or -328 ... +2372°F |
| <b>Type S (Pt10Rh-Pt):</b>          | -50.0 ... +999.9°C or -50 ... +1768°C   | -58.0 ... +999.9°F or -58 ... +3214°F   |
| <b>Type T (Cu-CuNi):</b>            | -120.0 ... +400.0°C or -220 ... +400°C  | -184.0 ... +752.0°F or -364 ... +752°F  |
| <b>Accuracy:</b>                    | (for thermoelements acc. to DIN EN 60584) ±1 digit (at nominal temperature)   |   |
| <b>Type K: -199.9 ... +999.9°C:</b> | ±0.03% of m.v. ±0.05%FS (T≥-60°C); ±0.2% of m.v. ±0.05%FS (T<-60°C)   |   |
| -220 ... +1372°C:                   | ±0.08% of m.v. ±0.1%FS (T≥-100°C); ±1°C ±0.1%FS (T<-100°C)  |   |
| <b>Type J: -120.0 ... +700.0°C:</b> | ±0.03% of m.v. ±0.08%FS (T≥-80°C); ±0.2% of m.v. ±0.08%FS (T<-80°C)   |   |
| -200 ... +1100°C:                   | ±0.08% of m.v. ±0.1%FS (T≥-150°C); ±1°C ±0.1%FS (T<-150°C)  |   |
| <b>Type N: -199.9 ... +999.9°C:</b> | ±0.03% of m.v. ±0.05%FS (T≥-60°C); ±0.2% of m.v. ±0.05%FS (T<-60°C)   |   |
| -200 ... +1300°C:                   | ±0.08% of m.v. ±0.1%FS (T≥-100°C); ±1°C ±0.1%FS (T<-100°C)  |   |
| <b>Type S: 0.0 ... +999.9°C:</b>    | ±0.05% of m.v. ±0.08%FS (T≥200°C); ±0.2% of m.v. ±0.08%FS (T<200°C)   |   |
| -50 ... +1768°C:                    | ±0.1% of m.v. ±0.1%FS (T≥200°C); ±1°C ±0.1%FS (T<200°C)   |   |
| <b>Type T: -120.0 ... +400.0°C:</b> | ±0.03% of m.v. ±0.1%FS (T≥-70°C); ±0.2% of m.v. ±0.1%FS (T<-70°C)   |   |
| -220 ... +400°C:                    | ±1°C (T≥-100°C); ±1°C ±1 digit (T<-100°C)   |   |
| <b>Temperature drift:</b>           | 0.01%/K   |   |
| <b>Point of comparison:</b>         | ±0.3°C  |   |
| <b>Nominal temperature:</b>         | 25°C  |   |
| <b>Working temperature:</b>         | -25 to +50°C  |   |
| <b>Relative humidity:</b>           | 0 to 95 %RH (non-condensing)  |   |
| <b>Storage temperature:</b>         | -25 to +70°C  |   |
| <b>Sensor connection:</b>           | 2 connection pins for miniature DIN-plug type K   |   |
| <b>Display :</b>                    | 2 four digit LCDs (12.4mm high and 7 mm high) for temperature, and for min./ max values, hold function, etc. as well as additional pointing arrows.   |   |
| <b>Pushbuttons:</b>                 | 6 membrane keys altogether for on/off switch, selection of thermoelements, min. and max. value memory, hold-function etc.   |   |
| <b>Interface:</b>                   | serial interface (3.5mm jack), serial interface can be directly connected to RS232 interface of a PC  |   |
| <b>Power supply:</b>                | 9V-battery, type IEC 6F22 (included) or additional d.c. connector (internal pin Ø 1.9 mm) for external 10.5-12V direct voltage supply. <br>(suitable power supply: PS 100) |   |
| <b>Power consumption:</b>           | approx. 1.6 mA  |   |
| <b>Automatic-off-function:</b>      | Device will be automatically switched off if no key is pressed/no interface communication takes place for the time of the power-off delay. The power-off delay can be set to values between 1 and 120 min.; it can be completely deactivated.               |   |
| <b>Min-/max-value memory:</b>       | Both the max. and the min. value will be memorized for sensor 1, sensor 2 and the difference.   |   |
| <b>Hold-function:</b>               | Press button to store current values of sensor 1, sensor 2 and their difference   |   |
| <b>Min./max alarm:</b>              | min. and max. values set for measuring values of sensor 1, sensor 2, sensor 1 and 2 or temperature difference are constantly monitored  |   |
| <b>Housing dimensions:</b>          | 142 x 71 x 26 mm (L x W x D)<br>impact-resistant ABS plastic housing, membrane keyboard, transparent panel. Front side IP65, integrated pop-up clip for table top or suspended use.   |   |
| <b>Weight:</b>                      | approx. 155 g   |   |
| <b>EMC:</b>                         | The device corresponds to the essential protection ratings established in the Regulations of the Council for the Approximation of Legislation for the member countries regarding electromagnetic compatibility (2004/108/EG)<br>Additional fault: <1%       |   |

## 12 Technical service

### Web Resources

Visit the VWR's website at [www.vwr.com](http://www.vwr.com) for:

- Complete technical service contact information
- Access to VWR's Online Catalogue, and information about accessories and related products
- Additional product information and special offers

**Contact us** For information or technical assistance contact your local VWR representative or visit [www.vwr.com](http://www.vwr.com).

## 13 Warranty

VWR International warrants that this product will be free from defects in material and workmanship for a period of two (2) years from date of purchase. If a defect is present, VWR will, at its option, repair, replace, or refund the purchase price of this product at no charge to you, provided it is returned during the warranty period. This warranty does not apply if the product has been damaged by accident, abuse, misuse, or misapplication, or from ordinary wear and tear.

For your protection, items being returned must be insured against possible damage or loss. This warranty shall be limited to the replacement of defective products. IT IS EXPRESSLY AGREED THAT THIS WARRANTY WILL BE IN LIEU OF ALL WARRANTIES OF FITNESS AND IN LIEU OF THE WARRANTY OF MERCHANTABILITY.

## 14 Equipment disposal



This equipment is marked with the crossed out wheeled bin symbol to indicate that this equipment must not be disposed of with unsorted waste.

Instead it's your responsibility to correctly dispose of your equipment at lifecycle-end by handling it over to an authorized facility for separate collection and recycling. It's also your responsibility to decontaminate the equipment in case of biological, chemical and/or radiological contamination, so as to protect from health hazards the persons involved in the disposal and recycling of the equipment.

For more information about where you can drop off your waste of equipment, please contact your local dealer from whom you originally purchased this equipment.

By doing so, you will help to conserve natural and environmental resources and you will ensure that your equipment is recycled in a manner that protects human health.

Thank you

**KALIBRIERZERTIFIKAT  
CALIBRATION CERTIFICATE  
CERTIFICAT DE CALIBRATION**



**Gerät / Device / Modèle**

TD 241

**Seriennummer  
Serial Number  
N° Série**

00774

**Messgröße**

**Meas. Unit**

**Unité de mesure**

**Temperatur**

**temperature**

**température**

**Sollwert  
reference  
référence**

0,0 °C

**Istwert  
actual value  
actuel**

0,2 °C

**Toleranz  
tolerance  
tolérance**

± 0,4 °C

**Gerät optisch  
Visual inspection  
inspection visuelle**

OK



**Datum / Date / Date**

22.03.2016

**Prüfer / Inspector / Verificateur**

JZ

VWR International Europe BVBA  
Researchpark Haasrode 2020  
Geldenaaksebaan 464  
B-3001 Leuven